

Solvent Waste Containment







Have you ever been written up for improper collection of solvent waste?

Consider this: For the cost of an HPLC column, you could be compliant. Plus, not collecting solvent waste properly poses a risk to the people and property in your lab.

The Problem The Solution **Ensures compliance with** Does not comply with safety regulations for safety standards your laboratory no more fines or write ups No tubing connections into waste About the price of a standard HPLC Column X Parafilm ≠ vapor control Unstable bottle Vapor protections Safeguards against accidental spills Stable, Compliant Unstable 4L **Chrom Tech Solvent Amber Bottle Waste Container** System Kit



Choosing the right system for solvent waste doesn't have to be complicated.

How many HPLCs are in your lab and are they adjacent to each other? We can accommodate multiple HPLCs into one 6-port manifold.

■ TECH TIP

Polypropylene or Stainless Steel? Choose Polypropylene (EPDM O-Ring) quick connects for typical reversed phase applications (economical and recommended option). Choose Stainless Steel (PTFE O-Ring) quick connects for normal phase or GPC applications.

SAFETY WASTE CONTAINER KIT-2 OR 5 GALLON KITS

6-PORT MANIFOLD, CONNECTIONS, AND ACCESSORIES INCLUDED IN KIT

- Organized laboratory space
- Remove harmful gases protect lab personnel from adverse health effects
- Tubing connections for the following tubing sizes:

1/16" OD (Qty 2) 1/8" OD (Qty 2) 1/4" ID (Qty 2) Plugs to close off any unused ports (Qty 4)










CHROM TECH SOLVENT WASTE CONTAINER COMPLETE SYSTEM KIT			
PART NO DESCRIPTION			
SAFETY-2PP-E	2 gal PE Solvent Waste Containment System w/PP quick connects		
SAFETY-5PP-E	SAFETY-5PP-E 5 gal PE Solvent Waste Containment System w/PP quick connects		
SAFETY-2SS-E	2 gal PE Solvent Waste Containment System w/SS quick connects		
SAFETY-5SS-E	5 gal PE Solvent Waste Containment System w/SS quick connects		


SAFETY KIT ACCESSORIES (NOT INCLUDED IN KIT)

IDEX BOTTLE CAPS, PLUGS & ADAPTERS

- Safely store your HPLC solvents
- Ports for 1/8" tubing and a luer taper
- Plugs and adapters for unused ports

IDEX BOTTLE CAPS					
	PORT				
PART NO	QTY	PORT STYLE	MATERIALS		
IDEX GL38 I	IDEX GL38 BOTTLE CAPS				
A-622	3	(2) 1/8" OD Tubing, (1) Luer Taper	PTFE, Polypropylene		
IDEX GL45 I	IDEX GL45 BOTTLE CAPS				
A-620	3	(2) 1/8" OD Tubing , (1) Luer Taper	PTFE, Polypropylene		
SAFETY A	SAFETY ACCESSORIES				
PART NO	DESCF	RIPTION			
JT-84003	HPLC	Can Spill Basin with 5 g spill capacity	/		
.IT-08500	Antistatic flexible wire with dual alligator clips 3'				

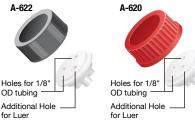

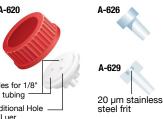


IDEX PLUGS & ADAPTERS		
PART NO	DESCRIPTION	
A-626	Bottle Cap Plug for luer hole, UHMWPE	
A-628	Bottle Cap Plug for 1/16", 1/8" or 3/16" hole, UHMWPE	
A-629	Filter Bottle Cap Plug for 1/16", 1/8" or 3/16" hole, UHMWPE with 20 μm stainless steel frit	





Do you need a better way to empty your safety disposal cans?

We make waste disposal SAFE and EASY with the All-Flo Pump available now through Chrom Tech!

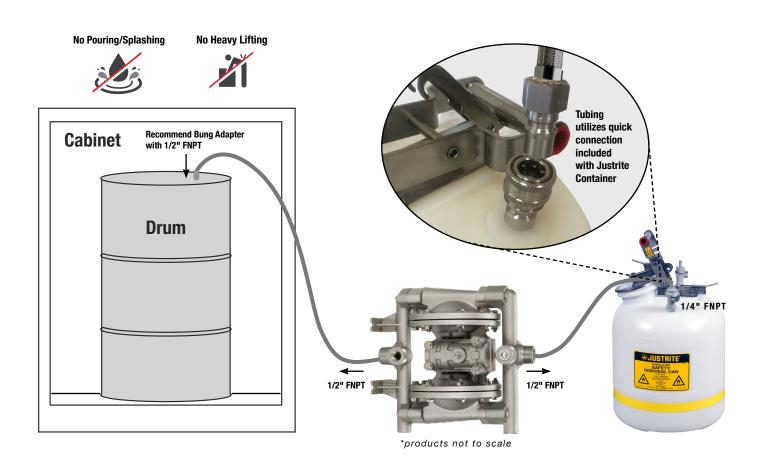

Tubing

Static Conductive Tubing PTFE-Lined, SS braided

Pump

A050-NP3-TT3T-S70 -All-Flo pump SS/PTFE 1/2" FNPT threads **Safety Can**

JT-12752-CC1 - 2 gal PP solvent waste safety can





SAFETY WASTE CONTAINERS: 2- OR 5-GALLON KITS

- · All use SS solvent pick up adapter
- · Available with PP or SS fitting adapters

ORGANIC VAPOR FILTER

- · Reusable adapter, disposable filters
- Remove harmful gases protect lab personnel from adverse health effects

6 PORT WASTE MANIFOLD WITH QUICK CONNECTIONS

- For various tubing sizes
- · Organized laboratory space



CHROM TECH SOLVENT WASTE SYSTEMS					
POLYPROPYLENE STAINLESS STEEL		DESCRIPTION & CAPACITY	DIMENSIONS	TUBING*	
JT-12752-CC1	JT-12752-CC3	2 gal (7.5L) Polyethylene safety disposal can w/SS fitting suction option assy	12" OD x 14.75" H	3/8"	
JT-12755-CC1	JT-12755-CC3	5 gal (19L) Polyethylene safety disposal can w/SS fitting suction option assy	12" OD x 20" H	3/8"	
MA-601	MA-603	6-Port Polypropylene waste manifold (5) 1/4" NPT, (1) 1/2" NPT			
FA-017	FA-019	Justrite can filter adapter elbow (to be used with CT-SCF-1)			
UNIV	ERSAL	DESCRIPTION & CAPACITY			
CT-S	SCF-5	Organic vapor filter, 5/pk			
EF	-102	Color Indicator for organic vapor filters, ea			
AB	-114	Barbed tubing connection, 1/4" ID, 1/4" NPT, PP			
AC	-150	Tubing connection, 1/16" OD, 1/4" NPT PP			
AC-155		Tubing connection, 1/8" OD, 1/4" NPT PP			
PL-101		Plug, 1/4" NPT, PP			
PL-102		Plug, 1/2" NPT, PP			

ALL-FLO STAINLESS STEEL VACUUM PUMP - A050

- Air-operated diaphragm pump—no electricity, eliminating the risk of sparks
- · Wetted materials: Stainless Steel and PTFE



■ TECH TIP

List of Wetted Components:PTFE Diaphragm, PTFE Ball, SS Seat, PTFE O-Ring, SS Hardware

ALL-FLO STAINLESS STEEL VACUUM PUMP- A050			
PART NO	DESCRIPTION		
A050-NP3-TT3T-S70	ALL-FLO pump, 1/2" female NPT ports, SS & PTFE components (see Tech Tip)		
AWE-050-TT3T	ALL-FLO pump maintenance kit		
SS-TC8PM4PM8-72	C8PM4PM8-72 Static conductive tubing 6' length, PTFE-Lined SS braided tube with 1/4" Male NPT to 1/2" Male NPT terminations (waste container to All-Flo pump)		
SS-TC8PM8PM8-72 Static conductive tubing 6' length, PTFE-Lined SS braided tube with ½" Male NPT to ½" Male NPT terminations (All-Flo pump to drum*)			

^{*}Need 1/2" Female NPT bung drum adapter



Disposal Cans (No Kit)

■ TECH TIP

Polypropylene quick-disconnect fittings are typically used with reverse phased solvents i.e. ACN, Methanol.

Stainless Steel/PTFE quick-disconnect fittings must be used with normal phase and GPC solvents i.e. Hexane, Heptane and THF.





Suction Option Assembly: Internal 1/4" (6 mm) ID polytetrafluoroethylene (PTFE) tubing with stainless steel fitting is installed below disconnect (of your specification) to allow for safe dispensing of fluid out of the can. NOTE: must be used under suction only.


FA-017

REPLACEMENT PARTS & ACCESSORIES FOR CHROM TECH SOLVENT WASTE CONTAINER SYSTEM KIT					
SAFETY DISPOSAL CANS					
DESCRIPTION & CAPACITY	DIMENSIONS	FITTINGS	0-RING	TUBING*	
1 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring	5.25" W x 12.75" H x 7.75" D	PP	EPDM	3/8"	
2 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring	12" OD x 14.75" H	PP	EPDM	3/8"	
5 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring	12" OD x 20" H	PP	EPDM	3/8"	
2 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring	12" OD x 14.75" H	SS	PTFE	3/8"	
5 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring	12" OD x 20" H	SS	PTFE	3/8"	
2 gal (7.5L) Polyethylene Safety Disposal Can w/SS fitting suction option assy	12" OD x 14.75" H	PP	EPDM	3/8"	
5 gal (19L) Polyethylene Safety Disposal Can w/SS fitting suction option assy	12" OD x 20" H	PP	EPDM	3/8"	
	DESCRIPTION & CAPACITY 1 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 2 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 5 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 2 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 5 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 2 gal (7.5L) Polyethylene Safety Disposal Can w/SS fitting suction option assy	DIMENSIONS 1 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 2 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 5.25" W x 12.75" H x 7.75" D 2 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 12" OD x 14.75" H 5 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 12" OD x 20" H 2 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 12" OD x 14.75" H 5 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 12" OD x 20" H 2 gal (7.5L) Polyethylene Safety Disposal Can w/SS fitting suction option assy 12" OD x 14.75" H	DIMENSIONS FITTINGS 1 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 5.25" W x 12.75" H x 7.75" D PP 2 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 12" OD x 14.75" H PP 5 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 12" OD x 20" H PP 2 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 12" OD x 14.75" H SS 5 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 12" OD x 20" H SS 2 gal (7.5L) Polyethylene Safety Disposal Can w/SS fitting suction option assy 12" OD x 14.75" H PP	DIMENSIONS FITTINGS O-RING 1 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 5.25" W x 12.75" H x 7.75" D PP EPDM 2 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 12" OD x 14.75" H PP EPDM 5 gal Polyethylene Safety Disposal Can, PP fittings, EPDM O-ring 12" OD x 20" H PP EPDM 2 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 12" OD x 14.75" H SS PTFE 5 gal Polyethylene Safety Disposal Can, SS fittings, PTFE O-ring 12" OD x 20" H SS PTFE 2 gal (7.5L) Polyethylene Safety Disposal Can w/SS fitting suction option assy 12" OD x 14.75" H PP EPDM	

^{*}Tubing is not supplied with individual containers

Chemical Emissions Filters & Adapters

- Minimize odors by adsorption before they enter the lab environment
- Constructed of activated coconut shell carbon and silica desiccant, with Polypropylene housing

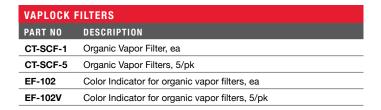



EF-102

JUSTRITE CAN ADAPTERS		
PART NO	DESCRIPTION	
FA-010	Justrite Can Adapter, EPDM, PP	
FA-015	Justrite Can Adapter, SS, PTFE	
FA-017	Justrite Can Adapter Elbow, EPDM, PP	
FA-019	Justrite Can Adapter Elbow, SS, PTFE	



Popular Fittings & Tubing



POPULAR MAN	NIFOLD FITTINGS & TUBING				
PART NO	TUBING	THREAD	MATERIAL	COLOR	QTY
HARDWALL TUB	ING ADAPTERS				
AC-150	1/16" OD	1/4" NPT	Polypropylene, ETFE	Green	ea
XP-201	1/16" OD	1/4-28	Delrin, ETFE	Black, Blue	ea
AC-155	1/8" OD	1/4" NPT	Polypropylene, ETFE	Blue	ea
XP-301	1/8" OD	1/4-28	Delrin, ETFE	Black, Yellow	ea
BARBED SOFTW	ALL TUBING ADAPTERS				
AB-114X	1/4" ID	1/4" NPT	Polypropylene		10/pk
AB-204X	1/4" ID	1/2" NPT	Polypropylene		10/pk
SOFTWALL ADA	PTER PLUGS FOR 1/4" AND 1/2" NP	T PORTS			
PL-101X	Plug	1/4" NPT	Polypropylene		10/pk
PL-102X	Plug	1/2" NPT	Polypropylene		10/pk
TUBING *Lengths	s greater than 10' available. Specify ler	ngth at time of order.			
T-10202	1/16" OD, 10'* length	-	FEP	Natural	ea
T-10201	1/8" OD, 10'* length	-	FEP	Natural	ea

HPLC Waste Containment Accessories

HPLC WAST	E CONTAINMENT ACCESSORIES
PART NO	DESCRIPTION
SF-003	Safety Waste Funnel, 4" diameter, PP disconnect
SF-005	Safety Waste Funnel, 4" diameter, SS disconnect
JT-84003	Safety Spill Basin for HPLC waste container
JT-06500	Antistatic Wire, dual alligator clips
JT-06530	Fill Pop Up Gauge







The Largest Chromatography Portfolio in the Industry



















































Chrom Tech, Inc.

5995 W 149th St Ste 102 Apple Valley, MN 55124 **TEL** 952-431-6000 sales@chromtech.com
online ordering | www.chromtech.com

CT-Lit 2302 rev1